ISCO 65X Syringe Pump



Brand: Teledyne ISCO Product Code: isco-65x Availability: Call for availability

Description

The 65x syringe pump provides precise, predictable flow and pressure control at flow rates from sub-microliter to 25 mL/min. Exceptional low-flow stability at up to 20,000 psi (1,390 bar) makes this the ideal pump for high pressure applications from deep well core analysis and polymer extrusions. The 65x pumps to the lowest flow rates with zero pulsation available on the market! The SyriXus controller has backlighted keypad and 4-line LCD display, as well as built-in and optional interfaces for computer control and other devices. Programming is easy and flexible, with instant access to menu screens even when the pump is running. This allows you to change operating parameters on the fly. Control the pump remotely using driver for LabVIEW[™] software, in addition to pump controller. The 65x pump module has a single-ratio drive train with autolubricating gears for long life and low maintenance. Special fittings are used to prevent leaks and ensure safety at maximum pressure. Wetted materials are compatible with most aqueous and organic liquids, corrosive solutions, heated fluids, liquefied gases. Continuous pumping is available in either single or dual pumping system controlled via air valves, offering zero pulsation for unlimited delivery. SyriXus pumps for proven reliable design. Known for years of performance, it can delivery fluid or pressurized gas for long term experiments. Or hold pressure constant in the system at a simple click of the button. Features</h2> 65x pump has 1/4" F-250 fittings for secure highpressure connections. Rugged single-ratio drive train has auto-lubricating gears to assure long life and low maintenance. SyriXus syringe pump is bench tested at the factory.

Key" controller. All ISCO SyriXus pumps use the same controller that can operate any four pump modules, either independently or together. Operating modes include: Single pump independent constant flow or pressure Single pump flow - or pressure - ramp programming dispensing Dual pump continuous constant flow or pressure >Dual pump concentration gradients>Four-pump configured systems

RS-232/485 serial interface for computer control or monitoring of operating parameters. SyriXus pumps have mechanical and electrical fail-safes, in addition to the factory default or user-set limits. standard material for cylinders, pistons, and caps - is an austenitic stainless steel with outstanding strength and corrosion resistance. heavy-duty, graphite-impregnated Teflon, providing very long life under harsh operating conditions. experiments requiring precision control nanobore capillary packed columns recovery studies in deep-well reservoirs chemical process development, catalyst evaluation, plastic formation <h2>Options and Accessories</h2> Temperature control jacket (cylinder cooling for filling with liquefied gas) programming software from National Instruments rate and pressure control.?? Capacity: 67.97 mL ? Flow Range: 0.00001 to 25 mL/min Flow Accuracy: + 0.3% of setpointPisplacement Resolution: 2.55 nL/step ?Press Range: 10 to 20,000 psi (?0.7 to 1,390 bar) Standard Pressure Accuracy: 0.5% Full ScaleOptional Pressure Accuracy: 0.1% Full Scale ?Wetted Materials: Nitronic 50, PTFE, Hastelloy C-276 ?Plumbing Ports: 1/4", F250C ?Operating Temperature: 5 to 40 °C Ambient VAC, 234 VAC, 50/60 Hz ?Dimensions (H x W x D): 39.8 x 10.7 x 18.4 inch (101 x 27 x 47 cm) ?Weight: ?Pump - 72.3 lb (32.8 kg), Controller -6.5 lb (3.0 kg?)